

CELLOTION®

Cell wash and recovery solution

- Serum-Free
- Chemically Defined

Cat#: 11898

Qty: 100ml


Expiry Date: 3 years from manufacturing date (see label)



CELLOTION® is a substitution for general-use cell wash solution such as PBS buffer, saline or culture medium with standards cell wash procedures.

Example of use: Monocytes isolation with CELLOTION® Cell Wash and Recovery Solution.

1. Gently dilute the peripheral blood or umbilical cord blood (UCB) with CELLOTION® Cell Wash and Recovery Solution 1:1.
2. Gently add Ficoll-hypaque solution to the mixture 1:2. Do not mix and the blood should remain on top.)
3. Centrifuge the tube for 30minutes at 1,800rpm, at room temperature.
4. Collect the buffy coat (white blood cells ring) carefully with a sterile pipette tip without touching the Ficoll-hypaque solution.
5. Gently dilute the cells with CELLOTION®, then wash for 2 - 3 times in CELLOTION®.
6. Count the number of cells.

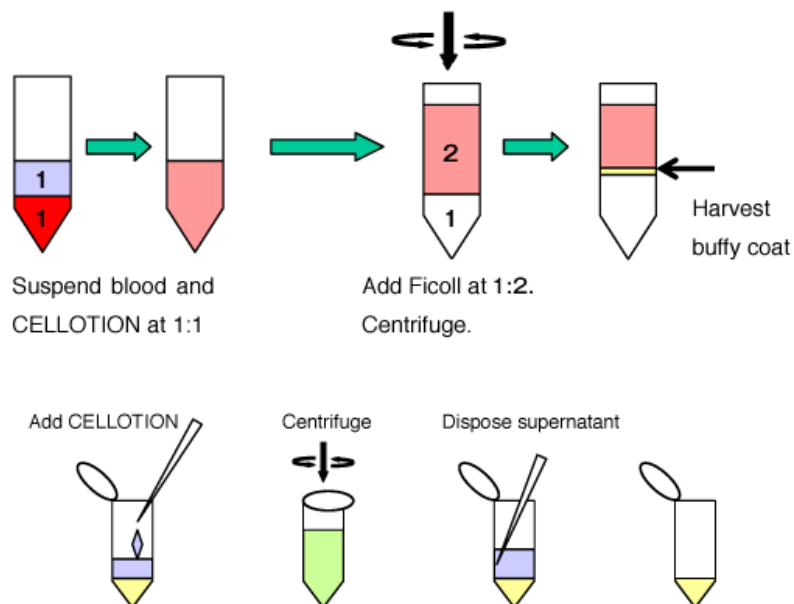

AMS BIO | www.amsbio.com | info@amsbio.com

 **UK & Rest of the World**
184 Park Drive, Milton Park
Abingdon OX14 4SE, UK
T: +44 (0)1235 828 200
F: +44 (0) 1235 820 482

 **North America**
1035 Cambridge Street,
Cambridge, MA 02141
T: +1 (617) 945-5033 or
T: +1 (800) 987-0985
F: +1 (617) 945-8218

 **Germany**
Bockenheimer Landstr. 17/19
60325 Frankfurt/Main
T: +49 (0) 69 779099
F: +49 (0) 69 13376880

 **Switzerland**
Centro Nord-Sud 2E
CH-6934 Bioggio-Lugano
T: +41(0) 91 604 55 22
F: +41(0) 91 605 17 85

Product Range:

Description	Pack Size
CELLBANKER 1 - Serum Containing	20 ml
CELLBANKER 1 - Serum Containing	100 ml
CELLBANKER 2 - Serum Free	20 ml
CELLBANKER 2 - Serum Free	100 ml
STEM-CELLBANKER® - GMP	20 ml
STEM-CELLBANKER® - GMP	100 ml
STEM-CELLBANKER® - GMP - DMSO Free	20 ml
STEM-CELLBANKER® - GMP - DMSO Free	100 ml
CELLOTION cell wash solution	100 ml

For research use only; not for use as a diagnostic.

AMSBIO | www.amsbio.com | info@amsbio.com

 **UK & Rest of the World**
184 Park Drive, Milton Park
Abingdon OX14 4SE, UK
T: +44 (0)1235 828 200
F: +44 (0) 1235 820 482

 **North America**
1035 Cambridge Street,
Cambridge, MA 02141
T: +1 (617) 945-5033 or
T: +1 (800) 987-0985
F: +1 (617) 945-8218

 **Germany**
Bockenheimer Landstr. 17/19
60325 Frankfurt/Main
T: +49 (0) 69 779099
F: +49 (0) 69 13376880

 **Switzerland**
Centro Nord-Sud 2E
CH-6934 Bioggio-Lugano
T: +41(0) 91 604 55 22
F: +41(0) 91 605 17 85