

Cell Culture Grade Bovine Type I Collagen

Catalog # 1202

For Research Use Only - Not Human or Therapeutic Use

Collagen, the most abundant protein in vertebrates, is observed in skin, cartilage, bone, intervertebral discs, blood vessels, tendons, ligaments, corneas, and is the main component of the extracellular matrix (ECM). It is identified at least 29 distinct collagen, type I collagen consists of two identical alpha 1 chains and one distinct alpha 2 chain (1), forming a triple helix formation known as tropocollagen. This triple-stranded helical conformation increases structural strength and resistance to enzymatic degradation, and plays a key role in assembly of the ECM.

Collagen is useful for facilitating tissue regeneration and/or site-specific drug delivery (2) because of its additional properties such as, low antigenicity, low toxicity, high water solubility, and high biodegradability. In particular, type I collagen binds integrins (3,4) facilitating cell migration (5), attachment (4,6), and proliferation and differentiation (3,6). Although atelocollagen, in which the telopeptides on the N- and C-terminals are removed from tropocollagen by pepsin digestion, is widely used in industrial purposes, tropocollagen may be beneficial as the native collagen scaffold containing cross-linked telopeptides.

We provide an acid soluble type I tropocollagen solution which can be used for traditional two-dimensional (2D) systems as well as a scaffold in three dimensional (3D) gels for simulating cell growth in fibroblasts (7,8) and chondrocytes (9). To determine the individual types of collagen, "tips for collagen solubilization protocols" are available, as well as ELISA kits for type I collagen and type II collagen. Please refer to www.amsbio.com or contact us for more details.

DESCRIPTION: Acid soluble bovine type I collagen solution

QUANTITY: 4 mg/ml x 12.5 ml, sterile-filtered

FORM: Dissolved in 0.01M HCl

PURITY: >95% type I collagen

SOURCE: Bovine

ENDOTOXIN: Less than 1 EU/ml

STORAGE TEMPERATURE: 4°C

STABILITY: 6 months

PROTOCOL: A) Plate Coating Procedures

Note: An optimal coating condition is required for your culture system.

1. Dilute the 4 mg/ml collagen with 0.02M HCl at 50 to 100 µg/ml.
2. Mix the diluted solution gently.
3. Add an appropriate volume of diluted collagen solution into wells or plates.

Note: Ensure the entire surface is coated.

4. Incubate at room temperature or 37°C for 1-2 hours.

AMSBIO | www.amsbio.com | info@amsbio.com



UK & Rest of the World
184 Park Drive, Milton Park
Abingdon OX14 4SE, UK
T: +44 (0)1235 828 200
F: +44 (0) 1235 820 482



North America
1035 Cambridge Street,
Cambridge, MA 02141
T: +1 (617) 945-5033 or
T: +1 (800) 987-0985
F: +1 (617) 945-8218



Germany
Bockenheimer Landstr. 17/19
60325 Frankfurt/Main
T: +49 (0) 69 779099
F: +49 (0) 69 13376880



Switzerland
Centro Nord-Sud 2E
CH-6934 Bioggio-Lugano
T: +41(0) 91 604 55 22
F: +41(0) 91 605 17 85

5. Remove all solution.
6. Rinse coated surfaces carefully with culture media or PBS.
7. The coated well or plates can be stored at 2-8°C or air dried if sterility is maintained.

B. Three-Dimensional (3D) Gel Preparation Procedures

1. Dilute the collagen solution with equal volume of cold sterilized PBS (final 2 mg/ml).
2. Add an appropriate volume of the diluted collagen solution in wells or plates.
3. Incubate at 37°C for 30-60 minutes.
4. The gel can be stored at 2-8°C or dried if sterility is maintained.

REFERENCES:

1. Shoulders, M. D. & Raines, R. T. Collagen structure and stability. Annual review of biochemistry 78, 929-958 (2009).
2. Friess, W. Collagen-biomaterial for drug delivery. European journal of pharmaceutics and biopharmaceutics: official journal of Arbeitsgemeinschaft fur Pharmazeutische Verfahrenstechnik e.V. 45, 113-136 (1998).
3. Taubenberger, A. V., Woodruff, M. A., Bai, H., Muller, D. J. & Hutmacher, D. W. The effect of unlocking RGD-motifs in collagen I on pre-osteoblast adhesion and differentiation. Biomaterials 31, 2827-2835 (2010).
4. Gullberg, D. et al. Analysis of alpha 1 beta 1, alpha 2 beta 1 and alpha 3 beta 1 integrins in cell-collagen interactions: identification of conformation dependent alpha 1 beta 1 binding sites in collagen type I. The EMBO journal 11, 3865-3873 (1992).
5. Bard, J. B. & Hay, E. D. The behavior of fibroblasts from the developing avian cornea. Morphology and movement in situ and in vitro. The Journal of cell biology 67, 400-418 (1975).
6. Kleinman, H. K., Klebe, R. J. & Martin, G. R. Role of collagenous matrices in the adhesion and growth of cells. The Journal of cell biology 88, 473-485 (1981).
7. Tomasek, J. J. & Hay, E. D. Analysis of the role of microfilaments and microtubules in acquisition of bipolarity and elongation of fibroblasts in hydrated collagen gels. The Journal of cell biology 99, 536-549 (1984).
8. Karamichos, D., Lakshman, N. & Petroll, W. M. Regulation of corneal fibroblast morphology and collagen reorganization by extracellular matrix mechanical properties. Investigative ophthalmology & visual science 48, 5030-5037 (2007).
9. Oliveira, S. M. et al. An improved collagen scaffold for skeletal regeneration. Journal of biomedical materials research. Part A 94, 371-379 (2010).

AMSBIO | www.amsbio.com | info@amsbio.com



UK & Rest of the World
184 Park Drive, Milton Park
Abingdon OX14 4SE, UK
T: +44 (0)1235 828 200
F: +44 (0) 1235 820 482



North America
1035 Cambridge Street,
Cambridge, MA 02141
T: +1 (617) 945-5033 or
T: +1 (800) 987-0985
F: +1 (617) 945-8218



Germany
Bockenheimer Landstr. 17/19
60325 Frankfurt/Main
T: +49 (0) 69 779099
F: +49 (0) 69 13376880



Switzerland
Centro Nord-Sud 2E
CH-6934 Bioggio-Lugano
T: +41(0) 91 604 55 22
F: +41(0) 91 605 17 85