

LS174T FFPE Cell Pellet

GENERAL INFORMATION

Product Name: LS174T FFPE Cell Pellet
 Reference Number: 3050-0410 Block
 3050-0420 Slide (4µm)
 3050-0430 FFPE scroll (20µm)
 Date of Manufacture: See product label
 Lot Number: See product label
 Intended Use: For research use only

*RECOMMENDED STABILITY RE-TESTING SCHEDULE:

Block	5 years
Scroll	1 year
Slide	1 year

*FFPE sample stability is not universally established and is biomarker and assay dependent. More frequent re-testing may be needed for certain labile biomolecules.

SAFETY AND PRECAUTIONS

This product does not contain hazardous material. Wear appropriate personal protective equipment (PPE) when handling reagents and biological specimens.

RECOMMENDED PROCEDURES

Staining using FFPE slides:

1. Bake slides at 60°C for 30-60 min.
2. Deparaffinize 3 times in Xylene or Xylene substitute for 5 min each time.
3. Rinse 2 times in 100% ethanol for 1 min each.
4. Rehydrate in ethanol series (95% 1 min, 70% 1 min, distilled H₂O 2 times for 1 min each).
5. Proceed to staining protocol.

Biomolecule extraction from FFPE scrolls:

1. Add 1ml Xylene or Xylene substitute to each tube containing FFPE scrolls and vortex for 30 sec.
2. Centrifuge at full speed for 2 min at room temperature. Remove supernatant without disturbing the pellet.
3. Repeat step 1 and 2 for a total of 3 times
4. Add 1ml 100% ethanol and mix by vortexing.
5. Centrifuge at full speed for 2 min at room temperature. Remove supernatant without disturbing the pellet.
6. Repeat step 4 and 5.
7. Carefully remove any residual ethanol in the tube without disturbing the pellet.
8. Open the tube and dry at room temperature or 37°C for 30min. Ensure that ethanol has completely evaporated.
9. Proceed to extraction protocol.

DESCRIPTION

Cell Line: LS174T
 Tissue of Origin: Colon or rectum
 Culturing Condition: DMEM supplemented with 10% FBS at 37°C with 5% CO₂
 Fixation Condition: 10% neutral buffered formalin (NBF) for 24 hours at 24-27°C
 Product Format:
 Block: Paraffin embedded block.
 Pellet diameter: 5mm
 Slide: One unstained section mounted on Superfrost™ Plus slide. Section thickness: 5µm
 FFPE Scroll: One FFPE section in DNase/RNase free tube. Section thickness: 20µm

SCHEMATICS

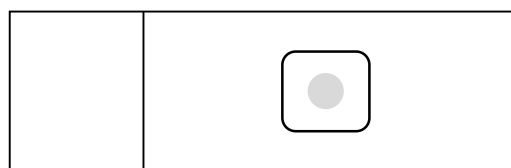


Illustration of an FFPE slide

STORAGE CONDITION

2-8 °C with desiccation (Recommended)
 For extended shelf life, store samples at -15 °C to -5 °C with desiccation.



AMSBIO | www.amsbio.com | info@amsbio.com

AMSBIO LLC
 USA & Canada 
 1035 Cambridge Street,
 Cambridge, MA 02141
 T: +1 (617) 945-5033 or
 T: +1 (800) 987-0985
 F: +1 (617) 945-8218

AMSBIO Europe BV
 EU 
 Berenkoog 41,
 1822 BH Alkmaar,
 Netherlands
 T: +31 (0) 72 8080244
 F: +31 (0) 72 8080142

AMS Biotechnology (Europe) Ltd
 UK & Rest of the World 
 184 Park Drive, Milton Park
 Abingdon OX14 4SE
 T: +44 (0) 1235 828 200
 F: +44 (0) 1235 820 482

AMS Biotechnology (Europe) Ltd
 Switzerland 
 Via Lisano 3,
 (CP.683)
 CH-6900 Massagno
 T: +41 (0) 91 604 55 22
 F: +41 (0) 91 605 17 85

AMSBIO Europe BV
 Deutschland 
 T: +49 (0) 69 779099
 F: +49 (0) 69 13376880