

## FITC-Dextran

Catalog # 4009

*For Research Use Only - Not Human or Therapeutic Use*

DESCRIPTION:	Fluorescein isothiocyanate labeled dextran
APPLICATIONS:	Use to assess the permeability of semi-permeable membranes either <i>in vivo</i> or <i>in vitro</i> (1, 2, 3).  Note: It is possible to simultaneously use FITC-dextran and TRITC-dextran of different sizes as FITC and TRITC fluoresce at different wavelengths.
QUANTITY:	FITC-dextran, 5 ml Note: 96-well plates (Catalog # 9045 are available upon request).
FORM:	25 mg/ml solution in PBS (green color).
MOLECULAR WEIGHT:	40 kDa
FLUORESCENT PROPERTIES:	Excitation: 490 nm, Emission: 520 nm
<i>IN VIVO</i> PROTOCOL:	<ol style="list-style-type: none"> <li>1. Fast mice 4 hours before the oral feeding, and for the duration of the experiment.</li> <li>2. Feed 20 ml/kg by oral gavage.</li> <li>3. Maintain fasting conditions and wait 3 hours (may vary depending on individual animals).</li> <li>4. Collect blood by retro-orbital bleeding, then spin and collect plasma. Dilute plasma 1:2 (or more) with PBS.</li> <li>5. Optional: Prepare a standard curve by making dilutions of the stock FITC-dextran with normal mouse plasma (12.5 ug/ml is a recommended starting point). <u>Caution: Protein in the samples may interfere and reduce the fluorescence intensity (FI).</u> Therefore, to ensure an accurate determination of FITC-dextran permeability, the Plasma to PBS ratio must be consistent throughout the standard curve, as well as with the dilution factor of the sample. Use normal mouse serum diluted in PBS for the blank.</li> <li>6. Transfer 50-100 ul of supernatant to a 96-well plate and read on a fluorescent plate reader.</li> </ol>
<i>IN VITRO</i> PROTOCOL:	Please refer to protocol of choice.
STORAGE TEMPERATURE:	4°C in the dark
STABILITY:	1 year

AMSBIO | [www.amsbio.com](http://www.amsbio.com) | [info@amsbio.com](mailto:info@amsbio.com)

**UK & Rest of the World**  
AMS Biotechnology (Europe) Ltd  
184 Park Drive, Milton Park  
Abingdon, UK  
T: +44 (0)1235 828 200  
F: +44 (0) 1235 820 482



**North America**  
amsbio LLC  
1035 Cambridge Street,  
Cambridge, MA 02141  
T: +1 617 945 5033 or  
T: +1 800 987 0985  
F: +1 617 945 8218



**Germany**  
AMS Biotechnology (Europe) Ltd  
Bockenheimer Landstr. 17/19  
60325 Frankfurt/Main  
T: +49 (0) 69 779099  
F: +49 (0) 69 13376880



**Switzerland**  
AMS Biotechnology (Europe) Ltd  
Centro Nord-Sud 2E  
CH-6934 Bioggio-Lugano  
T: +41(0) 91 604 55 22  
F: +41(0) 91 605 17 85

## REFERENCES:

1. M. Vijay-Kumar et al., Deletion of TLR5 results in spontaneous colitis in mice. J Clin Invest 117, 3909-3921 (2007).
2. Q. Wang et al., Intestinal permeability is reduced and IL-10 levels are increased in septic IL-6 knockout mice. Am J Physiol Regul Integr Comp Physiol 281, R1013-1023 (2001).
3. W. Huang et al., HMGB1 increases permeability of the endothelial cell monolayer via RAGE and Src family tyrosine kinase pathways. Inflammation 35, 350-362 (2012).

AMSBIO | [www.amsbio.com](http://www.amsbio.com) | [info@amsbio.com](mailto:info@amsbio.com)



**UK & Rest of the World**  
AMS Biotechnology (Europe) Ltd  
184 Park Drive, Milton Park  
Abingdon, UK  
T: +44 (0)1235 828 200  
F: +44 (0) 1235 820 482



**North America**  
amsbio LLC  
1035 Cambridge Street,  
Cambridge, MA 02141  
T: +1 617 945 5033 or  
T: +1 800 987 0985  
F: +1 617 945 8218



**Germany**  
AMS Biotechnology (Europe) Ltd  
Bockenheimer Landstr. 17/19  
60325 Frankfurt/Main  
T: +49 (0) 69 779099  
F: +49 (0) 69 13376880



**Switzerland**  
AMS Biotechnology (Europe) Ltd  
Centro Nord-Sud 2E  
CH-6934 Bioggio-Lugano  
T: +41(0) 91 604 55 22  
F: +41(0) 91 605 17 85