

## Data Sheet

### ***TNKS1 Histone Ribosylation Assay Kit (Biotin-labeled NAD<sup>+</sup>)*** Catalog # AMS.80579

**DESCRIPTION:** The TNKS1 Histone Ribosylation Assay Kit (Biotin-labeled NAD<sup>+</sup>) is designed to measure Tankyrase 1 (TNKS1) activity for screening and profiling applications. TNKS1 catalyzes the NAD-dependent addition of poly(ADP-ribose) to the substrate proteins. The TNKS1 assay kit comes in a convenient 384-well format, with purified TNKS1 enzyme, histone mixture, and PARP assay buffer for 384 enzyme reactions. The key to the TNKS1 Histone Ribosylation Assay (Biotin-labeled NAD<sup>+</sup>) is the biotinylated substrate. With this kit, only three simple steps are required for TNKS1 reactions. First, histone proteins are coated on a 384-well plate. Next, the biotinylated TNKS1 substrate is incubated with an assay buffer that contains the TNKS1 enzyme. Finally, the plate is treated with streptavidin-HRP followed by addition of the HRP substrate to produce chemiluminescence that can be measured using a chemiluminescence reader.

#### COMPONENTS:

Catalog #	Component	Amount	Storage
80504	TNKS1	2 x 5 µg	-80°C
52029	5x histone mixture	2 x 1 ml	-80°C
80601	10x Assay Mixture Containing Biotinylated Substrate	2 x 300 µl	-80°C
80602	10x PARP assay buffer	2 x 1 ml	-20°C
	Blocking buffer	2 x 25 ml	+4°C
	Streptavidin-HRP	2 x 100 µl	+4°C
	HRP chemiluminescent substrate A (translucent bottle)	2 x 6 ml	+4°C
	HRP chemiluminescent substrate B (brown bottle)	2 x 6 ml	+4°C
	384-well plate	1	Room Temperature

#### MATERIALS AND INSTRUMENTS REQUIRED BUT NOT SUPPLIED:

1x PBS buffer

PBST buffer (1x PBS, containing 0.05% Tween-20)

Luminometer or fluorescent microplate reader capable of reading chemiluminescence

Adjustable micropipettor and sterile tips

Rotating or rocker platform

RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

[AMS BIO | www.amsbio.com](http://www.amsbio.com) | [info@amsbio.com](mailto:info@amsbio.com)

 **UK & Rest of the World**  
184 Park Drive, Milton Park  
Abingdon OX14 4SE, UK  
T: +44 (0)1235 828 200  
F: +44 (0) 1235 820 482

 **North America**  
1035 Cambridge Street,  
Cambridge, MA 02141  
T: +1 (617) 945-5033 or  
T: +1 (800) 987-0985  
F: +1 (617) 945-8218

 **Germany**  
Bockenheimer Landstr. 17/19  
60325 Frankfurt/Main  
T: +49 (0) 69 779099  
F: +49 (0) 69 13376880

 **Switzerland**  
Centro Nord-Sud 2E  
CH-6934 Bioggio-Lugano  
T: +41(0) 91 604 55 22  
F: +41(0) 91 605 17 85

**APPLICATIONS:** Great for studying enzyme kinetics and screening small molecular inhibitors for drug discovery and HTS applications.

**STABILITY:** Up to 1 year when stored as recommended.

**REFERENCE:**

Brown, J.A., Marala, R.B. *J. Pharmacol. Toxicol. Methods* 2002 **47**:137-41.

**Assay Protocol:**

***All samples and controls should be tested in duplicate.***

Coating the plate with the histone mixture:

- 1) Dilute 5x histone mixture 1:5 in PBS.
- 2) Add 25  $\mu$ l diluted histone mixture to each well and incubate overnight at 4°C.
- 3) Wash the plate 3 times with 100  $\mu$ l PBST buffer.
- 4) Block the wells by adding 100  $\mu$ l of Blocking buffer to every well. Incubate for 60 minutes at room temperature.
- 5) Wash the plate 3 times with 100  $\mu$ l PBST buffer

(Alternatively, the plate can be coated for 90 minutes at 37°C followed by 60 minutes blocking at room temperature. All washing steps should be the same.)

**Step 1: Ribosylation reaction**

- 1) Prepare the master mixture: N wells x (1.25  $\mu$ l **10x PARP assay buffer** + 1.25  $\mu$ l **10x PARP assay mixture** + 10  $\mu$ l H<sub>2</sub>O)
- 2) Thaw **TNKS1 enzyme** on ice. Upon first thaw, briefly spin tube containing enzyme to recover full content of the tube. Aliquot **TNKS1 enzyme** into single use aliquots. Store remaining undiluted enzyme in aliquots at -80°C immediately. *Note: TNKS1 enzyme is very sensitive to freeze/thaw cycles. Do not re-use thawed aliquots or diluted enzyme.*
- 3) Dilute **TNKS1 enzyme** in **1X PARP assay buffer** at 1.25 – 2.5 ng/ $\mu$ l (12.5 - 25 ng/10  $\mu$ l). Keep diluted enzyme on ice until use. Discard any unused diluted enzyme after use.

Add 12.5  $\mu$ l of master mixture to each well designated for the “Positive Control”, “Test Inhibitor”, and “Blank”. For the “Substrate Control”, add 1.25  $\mu$ l **10x PARP assay buffer** + 11.25  $\mu$ l H<sub>2</sub>O.

RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.


	Blank	Positive Control	Substrate Control	Test Inhibitor
10X PARP Assay Buffer	1.25 $\mu$ l	1.25 $\mu$ l	1.25 $\mu$ l	1.25 $\mu$ l
10X assay mixture	1.25 $\mu$ l	1.25 $\mu$ l	-	1.25 $\mu$ l
H <sub>2</sub> O	10 $\mu$ l	10 $\mu$ l	11.25 $\mu$ l	10 $\mu$ l
Test Inhibitor	-	-	-	2.5 $\mu$ l
Inhibitor buffer (no inhibitor)	2.5 $\mu$ l	2.5 $\mu$ l	2.5 $\mu$ l	
1x PARP buffer	10 $\mu$ l	-	-	-
TNKS1 (~ 2 ng/ $\mu$ l)	-	10 $\mu$ l	10 $\mu$ l	10 $\mu$ l
<b>Total</b>	<b>25 <math>\mu</math>l</b>	<b>25 <math>\mu</math>l</b>	<b>25 <math>\mu</math>l</b>	<b>25 <math>\mu</math>l</b>

- 4) Add 2.5  $\mu$ l of inhibitor solution to each well designated "Test Inhibitor". For the "Positive Control", "Substrate Control", and "Blank", add 2.5  $\mu$ l of the same solution without inhibitor (inhibitor buffer).
- 5) Add 10  $\mu$ l of **1x PARP buffer** to the well designated "Blank".
- 6) Initiate the reactions by adding 10  $\mu$ l of diluted **TNKS1** prepared as described above. Incubate the reactions for 1 hour at room temperature.
- 7) Wash the plate 3 times with 100  $\mu$ l PBST per well.

### Step 3: Detection

- 1) Dilute **Streptavidin-HRP** 1:50 in Blocking buffer.
- 2) Add 25  $\mu$ l of diluted **Streptavidin-HRP** to each well. Incubate for 30 minutes at room temperature.
- 3) Wash three times with 100  $\mu$ l PBST buffer as above.
- 4) Just before use, mix on ice 25  $\mu$ l **HRP chemiluminescent substrate A** and 25  $\mu$ l **HRP chemiluminescent substrate B** and add 50  $\mu$ l per well. Discard any unused chemiluminescent reagent after use.
- 5) Incubate for 5 minutes at room temperature then read sample in a luminometer or microtiter-plate reader capable of reading chemiluminescence  
RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

**Example of Assay Results:**

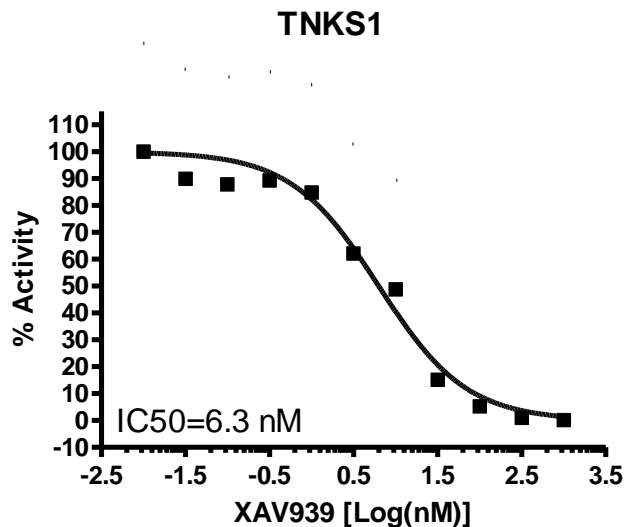

Inhibition of TNKS1 enzyme (Cat. # 80504) with XAV939 (Cat. # 27100), measured using the TNKS1 Histone Ribosylation Assay Kit (Biotin-labeled NAD<sup>+</sup>), AMSBIO (Cat. # 80579). Luminescence was measured using a Bio-Tek fluorescent microplate reader. *Data shown is lot-specific.*

**RELATED PRODUCTS:**

<u>Product</u>	<u>Cat. #</u>	<u>Size</u>
TNKS1 Histone Ribosylation Assay Kit (Antibody Detection)	80574	96 rxns.
TNKS2 Histone Ribosylation Assay Kit (Antibody Detection)	80576	96 rxns.
TNKS2 Histone Ribosylation Assay Kit (Biotin-labeled NAD <sup>+</sup> )	80578	96 rxns.
TNKS1 (PARP5A) Enzyme	80504	10 µg
TNKS2 (PARP5B), (667-end) Enzyme	80505	10 µg
TNKS2 (PARP5B), (849-end) Enzyme	80515	10 µg
PARP5b (TNKS2) Assay Kit	80576	96 rxns.
PARP6 Assay Kit	80556	32 rxns.
PARP1 Enzyme	80501	10 µg
PARP2 Enzyme	80502	10 µg
PARP3 Enzyme	80503	10 µg
PARP6 Enzyme	80506	10 µg

AMSBIO | [www.amsbio.com](http://www.amsbio.com) | [info@amsbio.com](mailto:info@amsbio.com)

 **UK & Rest of the World**  
 184 Park Drive, Milton Park  
 Abingdon OX14 4SE, UK  
 T: +44 (0)1235 828 200  
 F: +44 (0) 1235 820 482

 **North America**  
 1035 Cambridge Street,  
 Cambridge, MA 02141  
 T: +1 (617) 945-5033 or  
 T: +1 (800) 987-0985  
 F: +1 (617) 945-8218

 **Germany**  
 Bockenheimer Landstr. 17/19  
 60325 Frankfurt/Main  
 T: +49 (0) 69 779099  
 F: +49 (0) 69 13376880

 **Switzerland**  
 Centro Nord-Sud 2E  
 CH-6934 Bioggio-Lugano  
 T: +41(0) 91 604 55 22  
 F: +41(0) 91 605 17 85