

amsbio» Accelerate Your Discovery

Version 003

Store at 2-15°C

For research use only Not for use in diagnostic procedures

# iMatrix-221

Product No. AMS 892 061	$350~\mu{ m g}$
Product No. AMS.892 062	$1,050~\mu{ m g}$

Product description: iMatrix-221 is a recombinant human laminin-221 E8 fragment protein expressed in Chinese Hamster Ovary (CHO)-S cells. iMatrix-221 contains the integrin-binding site of the laminin-221 molecule. iMatrix-221 is a useful cell culture substrate for proliferation and differentiation of cardiomyocytes and skeletal muscule cells. iMatrix-221 is also useful for the culture of other cells adhering to laminin-221.

Content: Recombinant human laminin-221  $\mathbf{E8}$ fragment protein in PBS(-)

Concentration: 0.5 mg/mL

Amount: 175 µg / 0.35 mL / tube Product No. 892 061, 892 066 350 µg / 2 tubes Product No. 892 062, 892 065 1,050 µg / 6 tubes

Storage: Store at 2°C to 15°C, protect from light.

**Expiration date:** The shelf life is 2 years from the date of manufacture. The expiration date is printed on the outer carton.

Activity: The dissociation constant (Kd) for the binding with integrin a7X261 is 10 nM or less.

Methods of use: By the following method, iMatrix-221 can be coated onto a culture vessel. The optimum coating density may differ by cell-type, cell-line, medium selected, or purpose. Insufficient coating density may result in the detachment of cells and varied cell conditions while the excessive coating density may lead to difficulty in detaching cells for passage.

Determine the optimal coating density.  $0.5 \,\mu\text{g/cm}^2$  is a standard but test between 0.1 and 1.5 µg/cm<sup>2</sup>.

- 1) Dilute iMatrix-221 with PBS(-). Use the diluted iMatrix-221 immediately. To coat with 0.5 µg/cm<sup>2</sup> onto a 6-well plate with 9.6 cm<sup>2</sup>/well, dilute 9.6 µL of iMatrix-221 with 2 mL of PBS(-) per well.
- 2) Place the diluted iMatrix-221 into a culture vessel and incubate either at 37°C for 1 h, or at room temperature for 3 h, or at 4°C overnight.
- Aspirate the coating solution. Then, immediately 3) seed your cells. Do not allow the coated surface to dry.

\*If you face difficulties in detaching cells for passage, re-adjust the conditions (e.g., reduce the coating density).

### **References**:

Nishiuchi R. et al. (2006), Matrix Biol. 25 (3): 189-97 Taniguchi Y. et al. (2009), J. Biol. Chem. 284 (12): 7820-31

Israeli-Rosenberg S. et al. (2014), Circ. Res. 114 (3): 572-86

Ishii K. et al. (2018), Stem Cell Reports 10 (2): 562-82

Caution: For research use only. Not intended for human use. In the event of accidental ingestion or contact with the eyes, immediately wash the affected area and seek medical attention.

Product information: Current information including references and Q&A are available at https://www.amsbio.com/imatrix-recombinantlaminin-series/

## **Designed by:** MATRIXOME, Inc.

3-2 Yamadaoka, Suita, Osaka 565-0871, Japan Institute for Protein Research, Osaka University

## Manufactured by: Nippi, Incorporated

1-1-1 Senju Midori-cho, Adachi, Tokyo 120-8601, Japan

AMSBIO | www.amsbio.com | info@amsbio.com

#### MSBIO LLC USA & Canada

1035 Cambridge Street, Cambridge, MA 02141 T: +1 (617) 945-5033 or T: +1 (800) 987-0985 F: +1 (617) 945-8218

#### AMSBIO Europe BV ΕU 1

Berenkoog 41, 1822 BH Alkmaar, Netherlands T: +31 (0) 72 8080244 F: +31 (0) 72 8080142

# AMS Biotechnology (Europe) Ltd UK & Rest of the World

184 Park Drive, Milton Park Abingdon OX14 4SE T: +44 (0) 1235 828 200 F: +44 (0) 1235 820 482

MS Biotechnology (Europe) Ltd Switzerland + Via Lisano 3 Via Lisano 3, (CP.683) CH-6900 Massagno T: +41 (0) 91 604 55 22 F: +41 (0) 91 605 17 85

# MSBIO Europe BV Deutschland

T: +49 (0) 69 779099 F: +49 (0) 69 13376880