



# Mimetix<sup>®</sup> 12-well plate

AMS.TECL-003-1X AMS.TECL-003-8X



Scaffold discs in the wells are removable, held in place by retaining rings and can be taken out for imaging purposes.

## Product Description:

In our Mimetix 12-well plate (supplied with lid) scaffold discs are held in place by retaining rings. Discs can be transferred to a new plate or mounted on a glass slide for imaging purposes.

## General Features:

- True 3D environment rather than a roughened 2D surface
- Minimal protocol adaption required to switch from 2D to 3D
- Compatible with fluorescence microscopy
- Scaffolds can be coated with materials to facilitate cell adhesion in low serum conditions
- Protocols for cell seeding, retrieval, assays, and imaging are available

## Scaffold Specifications:

- Material: medical-grade poly-L-lactide (PLLA)
- Orientation: Random, non-woven
- Thickness: 50 µm
- Fiber diameter: 4 µm (15-30 µm pores)
- Overall porosity: app. 80%
- Non-biodegradable in *in vitro* applications

## Plate specifications:

- Corning 12-well plate: see [https://www.corning.com/catalog/cls/documents/drawings/LSR00002\\_Multiwell\\_TC\\_Plates\\_12\\_Well.pdf](https://www.corning.com/catalog/cls/documents/drawings/LSR00002_Multiwell_TC_Plates_12_Well.pdf)
- Supplied with lid in individually-sealed plastic wrapping
- Treated with gamma or e-beam irradiation
- Store at room temperature in the dark
- Manufactured in the United Kingdom

AMSBIO | [www.amsbio.com](http://www.amsbio.com) | [info@amsbio.com](mailto:info@amsbio.com)



**UK & Rest of the World**  
AMS Biotechnology (Europe) Ltd  
184 Park Drive, Milton Park  
Abingdon, UK  
T: +44 (0)1235 828 200  
F: +44 (0) 1235 820 482



**North America**  
amsbio LLC  
1035 Cambridge Street,  
Cambridge, MA 02141  
T: +1 (617) 945-5033 or  
T: +1 (800) 987-0985  
F: +1 (617) 945-8218



**Germany**  
AMS Biotechnology (Europe) Ltd  
Bockenheimer Landstr. 17/19  
60325 Frankfurt/Main  
T: +49 (0) 69 779099  
F: +49 (0) 69 13376880



**Switzerland**  
AMS Biotechnology (Europe) Ltd  
Centro Nord-Sud 2E  
CH-6934 Bioggio-Lugano  
T: +41(0) 91 604 55 22  
F: +41(0) 91 605 17 85



# Mimetix<sup>®</sup> 12-well plate

### Precondition

The Mimetix scaffold needs to be wetted with ethanol in order to allow a cell suspension to access the pores.

- Add 1 mL 20% ethanol per well.
- Allow ethanol to soak into the membrane for 5 min, then aspirate ethanol carefully without touching the scaffold.

### Wash

- Wash scaffold twice with PBS.
- Leave scaffold in cell culture medium until cell seeding.

### Seed

These seeding densities are general guidelines only.

- Add 50,000-100,000 cells suspended in 1 mL cell culture medium.

### Exchange medium

- For long-term experiments semi-exchange the cell culture medium every 3 days.

AMSBIO | [www.amsbio.com](http://www.amsbio.com) | [info@amsbio.com](mailto:info@amsbio.com)



**UK & Rest of the World**  
AMS Biotechnology (Europe) Ltd  
184 Park Drive, Milton Park  
Abingdon, UK  
T: +44 (0)1235 828 200  
F: +44 (0) 1235 820 482



**North America**  
amsbio LLC  
1035 Cambridge Street,  
Cambridge, MA 02141  
T: +1 (617) 945-5033 or  
T: +1 (800) 987-0985  
F: +1 (617) 945-8218



**Germany**  
AMS Biotechnology (Europe) Ltd  
Bockenheimer Landstr. 17/19  
60325 Frankfurt/Main  
T: +49 (0) 69 779099  
F: +49 (0) 69 13376880



**Switzerland**  
AMS Biotechnology (Europe) Ltd  
Centro Nord-Sud 2E  
CH-6934 Bioggio-Lugano  
T: +41(0) 91 604 55 22  
F: +41(0) 91 605 17 85