



Specification Sheet

Select-HA™ 1000kEF
Catalogue No. HYA-1000kEF-1

Product Description

Select-HA™ is a hyaluronic acid (HA) preparation of uniform and narrow size distribution prepared by *in vitro* synthesis using recombinant *Pasteurella multocida* hyaluronan synthase¹. Select-HA™ 1000kEF is defined as Select-HA polymer with a specific average weight-average molecular mass (Mw) within the range of 800 kDa to 1200 kDa and the polymer contains minimal level of endotoxin. (Although there is lot to lot variation, the sizes of a particular lot are generally within $\pm 5\%$). **Please refer to Certificate of Analysis for the specific average Mw of the lot you have received.**

Quantity

The tube contains 1 mg of Select-HA™ supplied as lyophilized sodium salts.

Recommended Reconstitution Procedure

Centrifuge the tube for a few seconds to collect the Select-HA™ solids in the bottom of the tube. Carefully open and add desirable amount of sterile water or proper buffer directly to the bottom of the tube. Allow two hours at 4°C for sample rehydration and then mix well before use.

Storage

Store at -20°C or below. To avoid frequent freeze-thaw, aliquoting is recommended. Avoid contamination with microbes or HA-degrading enzymes.

References

1. Jing, W. and DeAngelis, P.L. (2004) Synchronized Chemoenzymatic Synthesis of Monodisperse Hyaluronan Polymers. *J. Biol. Chem.* **279**(40), 42345-42349.

For research use only



UK & Rest of World

184 Milton Park, Abingdon
OX14 4SE, Oxon, UK
Tel: +44 (0) 1235 828 200
Fax: +44 (0) 1235 820 482

Switzerland

Centro Nord-Sud 2E
CH-6934 Bioggio-Lugano
Tel: +41 (0) 91 604 55 22
Fax: +41 (0) 91 605 17 85

Deutschland

Bockenheimer Landstr. 17/19
60325 Frankfurt/Main
Tel: +49 (0) 69 779099
Fax: +49 (0) 69 13376880

United States

23591 El Toro Rd, Suite #167
Lake Forest, CA 92630
Tel: +1 855 267 2464
Fax: +1 949 768 8613

amsbio
info@amsbio.com