

# 96-well Serum/Plasma Glycerol Kit

## Free Glycerol Detection

Cat# SGA-1

INSTRUCTION MANUAL ZBM0043.00

### STORAGE CONDITIONS

---

- Reagents & Buffers: 4°C
- Glycerol Standard: -20°C
- Blank assay plates (96-well): Room Temperature

For *in vitro* Use Only

### LIMITED PRODUCT WARRANTY

This warranty limits our liability to replacement of this product. No other warranties of any kind, expressed or implied, including without limitation, implied warranties of merchantability or fitness for a particular purpose, are provided by Zen-Bio, Inc. Zen-Bio, Inc. shall have no liability for any direct, indirect, consequential, or incidental damages arising out of the use, the results of use, or the inability to use this product.

### ORDERING INFORMATION AND TECHNICAL SERVICES

The product accompanying this document is intended for research use only and is not intended for diagnostic purposes or for use in humans.



#### UK & Rest of World

184 Milton Park, Abingdon  
OX14 4SE, Oxon, UK  
Tel: +44 (0) 1235 828 200  
Fax: +44 (0) 1235 820 482

#### Switzerland

Centro Nord-Sud 2E  
CH-6934 Bioggio-Lugano  
Tel: +41 (0) 91 604 55 22  
Fax: +41 (0) 91 605 17 85

#### Deutschland

Bockenheimer Landstr. 17/19  
60325 Frankfurt/Main  
Tel: +49 (0) 69 779099  
Fax: +49 (0) 69 13376880

#### United States

23591 El Toro Rd, Suite #167  
Lake Forest, CA 92630  
Tel: + 1 800 987 0985  
Fax: + 1 949 265 7703


# INTRODUCTION

This kit is designed to accurately determine the amount of glycerol present in blood serum or plasma of humans, mice, rats, and other animals in a 96-well format for increased throughput analysis. Blood can be collected in plain evacuated tubes or in the presence of common anti-coagulants: sodium citrate, ammonium oxalate and EDTA. **NOTE: Heparin or Heparinized tubes should not be used because this will generate inaccurate readings.** Serum should be separated from clotted blood by centrifugation as soon as possible and may be stored frozen (-20°C) prior to analysis.

## PRINCIPLE OF THE ASSAY

### Detection of Free Glycerol

Glycerol present in sample is phosphorylated by adenosine triphosphate (ATP) forming glycerol-1-phosphate (G-1-P) and adenosine-5'-diphosphate (ADP) in the reaction catalyzed by glycerol kinase. G-1-P is then oxidized by glycerol phosphate oxidase to dihydroxyacetone phosphate (DAP) and hydrogen peroxide (H<sub>2</sub>O<sub>2</sub>). A quinoneimine dye is produced by the peroxidase catalyzed coupling of 4-aminoantipyrine (4-AAP) and sodium N-ethyl-N-(3-sulfopropyl)m-anisidine (ESPA) with H<sub>2</sub>O<sub>2</sub>, which shows an absorbance maximum at 540nm. The increase in absorbance at 540nm is directly proportional to glycerol concentration of the sample.

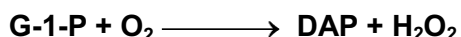

## ITEMS INCLUDED IN THE KIT

ITEM	DESCRIPTION	Cap Color	UNIT	QTY	STORAGE
Assay Plate, Plate A	96-well assay plate, blank	---	PLATE	2	-----
Dilution Buffer	12 ml	---	BOTTLE	1	4°C
Glycerol Reagent A	Reconstitute with 11.0 ml deionized water prior to use.	---	BOTTLE	1	4°C
Tray	For multi-channel pipettors, clear polyvinyl	CLEAR	EACH	2	-----
Glycerol standard	Glycerol @ 1mM [Dilute with 200 µl Dilution Buffer to make the 200 µM glycerol standard; see page 5 for recommended dilution scheme]	ORANGE	50 µl / VIAL	1	-20°C

### Other equipment/reagents required but not provided with the kit:

- Multi-channel Pipet , single channel pipet and pipet tips
- Plate reader with a filter of 540 nm
- Incubator at 37°C
- Large gauge needle
- Tubes for diluting standards



#### UK & Rest of World

184 Milton Park, Abingdon  
OX14 4SE, Oxon, UK  
Tel: +44 (0) 1235 828 200  
Fax: +44 (0) 1235 820 482

#### Switzerland

Centro Nord-Sud 2E  
CH-6934 Bioggio-Lugano  
Tel: +41 (0) 91 604 55 22  
Fax: +41 (0) 91 605 17 85

#### Deutschland

Bockenheimer Landstr. 17/19  
60325 Frankfurt/Main  
Tel: +49 (0) 69 779099  
Fax: +49 (0) 69 13376880

#### United States


23591 El Toro Rd, Suite #167  
Lake Forest, CA 92630  
Tel: + 1 800 987 0985  
Fax: + 1 949 265 7703

**amsbio**  
info@amsbio.com

# ASSAY PROCEDURE

## 1. Prepare the glycerol standards as follows:

Briefly spin down the contents of the glycerol standard tube before reconstitution. Pipette 200  $\mu\text{l}$  of Dilution Buffer into the 1 mM glycerol standard tube provided and mix well by vortexing. This produces a diluted stock glycerol standard of 200  $\mu\text{M}$ . Pipette 125  $\mu\text{l}$  of dilution buffer into 6 tubes (not provided). Using the newly diluted stock glycerol solution, prepare a dilution series as depicted below. Mix each new dilution thoroughly before proceeding to the next. The 200  $\mu\text{M}$  stock dilution serves as the highest standard, and the dilution buffer serves as the zero standard.

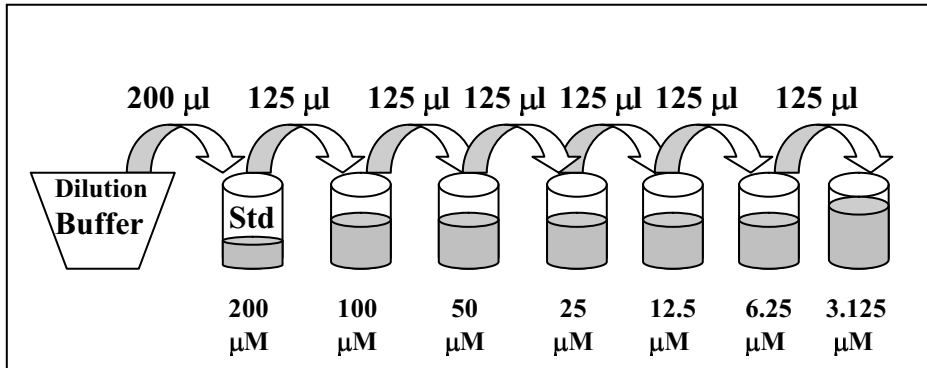

- Also at this time prepare the Glycerol Reagent A by adding 11.0 ml room temperature deionized water per bottle and gently invert. **DO NOT VORTEX!** Use a pipet to ensure that the powder is completely dissolved. Store at room temperature. If using a Reagent A solution previously prepared and stored at 2-8°C, also bring to room temperature. Make sure there is enough Reagent A from one solution to treat all the points in the assay. It may be necessary to combine solutions. Store in a light protected bottle. Reconstituted Glycerol Reagent A is stable for 60 days refrigerated (2-8°C); store any remaining solution refrigerated (2-8°C).
- Add 20  $\mu\text{l}$  (or 10 - 25  $\mu\text{l}$ ) of serum or plasma to a well of a blank plate for assessment of free glycerol. Add 30  $\mu\text{l}$  of dilution buffer to each well to total 50  $\mu\text{l}$  including serum or plasma sample. **THIS RESULTS IN A 2.5x DILUTION OF YOUR SAMPLE (20  $\mu\text{l}$  in 50  $\mu\text{l}$ )**. Add 50  $\mu\text{l}$  of each standard to empty wells (use another plate, if necessary).
- Add the reconstituted Glycerol Reagent A solution to one of the disposable trays provided in the kit. Add 50  $\mu\text{l}$  of Reagent A to each well of the glycerol plate. Gently, pipet up and down once to mix. Pop the bubbles using a large gauge needle or a clean pipet tip. The plate is then incubated at 25°C (room temperature) for 15 minutes.
- The optical density of each well is then measured at 540 nm.



### UK & Rest of World

184 Milton Park, Abingdon  
OX14 4SE, Oxon, UK  
Tel: +44 (0) 1235 828 200  
Fax: +44 (0) 1235 820 482

### Switzerland

Centro Nord-Sud 2E  
CH-6934 Bioggio-Lugano  
Tel: +41 (0) 91 604 55 22  
Fax: +41 (0) 91 605 17 85

### Deutschland

Bockenheimer Landstr. 17/19  
60325 Frankfurt/Main  
Tel: +49 (0) 69 779099  
Fax: +49 (0) 69 13376880

### United States

23591 El Toro Rd, Suite #167  
Lake Forest, CA 92630  
Tel: + 1 800 987 0985  
Fax: + 1 949 265 7703

**amsbio**  
info@amsbio.com

# GLYCEROL STANDARD CURVE


Generate standard curve: see example below

**DO NOT use this standard curve to generate your data. This is an example.]**

Subtract the OD value of the 0 $\mu$ M standard from all OD values including the standard curve.

Zero  
(blank) = .040

$\mu$ M Glycerol	OD	OD - blank
3.125	0.054	0.014
6.25	0.066	0.026
12.5	0.082	0.042
25	0.138	0.098
50	0.214	0.174
100	0.402	0.362
200	0.711	0.671

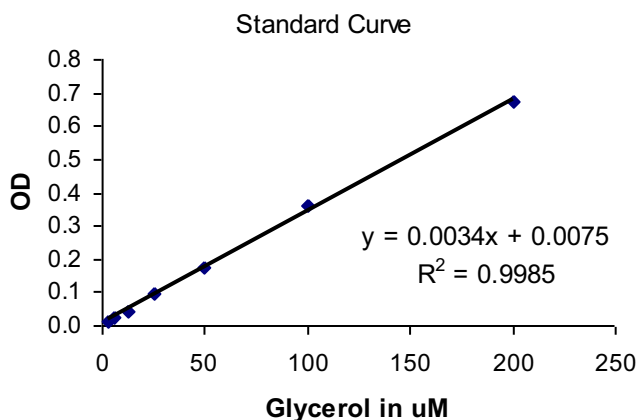

slope =	0.0034
intercept=	0.0075
$r^2$ =	0.9985

y = observed O.D. minus the blank

x = concentration of glycerol in  $\mu$ M

To calculate x for each y, (i.e. to change the observed O.D. into glycerol concentration) use the following equation:

y=(slope) times (x) plus intercept

y=mx+b so x=(y-b)/m

x=(y - 0.0075)/0.003 where 0.003= slope of the line and 0.0075= y intercept. Be careful to enter the proper sign for the y intercept value as it may be a negative number.

Any OD values greater than the highest standard (200  $\mu$ M), should be suspect. The compound should be re-assayed using a lower dose of the compound at treatment OR a dilute solution of the condition medium at the time of the assay.

**The  $R^2$  value should be equal or greater than 0.98 for the standard curve to be valid. Any  $R^2$  values below 0.98, must have the standard curve run again.**

Data are expressed as  $\mu$ M glycerol.

**REMEMBER TO ACCOUNT FOR THE DILUTION FACTOR IN STEP 3.**



UK & Rest of World

184 Milton Park, Abingdon  
OX14 4SE, Oxon, UK  
Tel: +44 (0) 1235 828 200  
Fax: +44 (0) 1235 820 482

Switzerland

Centro Nord-Sud 2E  
CH-6934 Bioggio-Lugano  
Tel: +41 (0) 91 604 55 22  
Fax: +41 (0) 91 605 17 85

Deutschland

Bockenheimer Landstr. 17/19  
60325 Frankfurt/Main  
Tel: +49 (0) 69 779099  
Fax: +49 (0) 69 13376880

United States

23591 El Toro Rd, Suite #167  
Lake Forest, CA 92630  
Tel: + 1 800 987 0985  
Fax: + 1 949 265 7703

**amsbio**  
info@amsbio.com

# APPENDIX A: PLATE LAYOUT

H	G	F	E	D	C	B	A	
								1
								2
								3
								4
								5
								6
								7
								8
								9
								10
								11
								12



**UK & Rest of World**  
 184 Milton Park, Abingdon  
 OX14 4SE, Oxon, UK  
 Tel: +44 (0) 1235 828 200  
 Fax: +44 (0) 1235 820 482

**Switzerland**  
 Centro Nord-Sud 2E  
 CH-6934 Bioggio-Lugano  
 Tel: +41 (0) 91 604 55 22  
 Fax: +41 (0) 91 605 17 85

**Deutschland**  
 Bockenheimer Landstr. 17/19  
 60325 Frankfurt/Main  
 Tel: +49 (0) 69 779099  
 Fax: +49 (0) 69 13376880

**United States**  
 23591 El Toro Rd, Suite #167  
 Lake Forest, CA 92630  
 Tel: + 1 800 987 0985  
 Fax: + 1 949 265 7703

**amsbio**  
 info@amsbio.com

# APPENDIX B: PROCEDURE FLOWCHART

## Glycerol Detection

Add 20  $\mu$ l/well test sample and 30  $\mu$ l/well dilution buffer to one of the blank assay plates provided.  
Add 50  $\mu$ l/well diluted standard curve to empty wells.



Reconstitute Glycerol Reagent A.  
Add 50  $\mu$ l/well.





Incubate 15 minutes @ room temperature.



Measure the optical density of each well at 540 nm using a spectrophotometer plate reader.

**Reminder: Sample was diluted in Step 3**

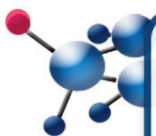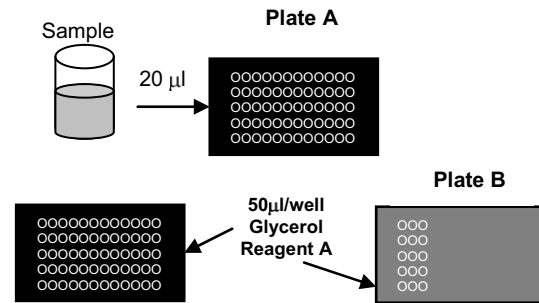

### UK & Rest of World

184 Milton Park, Abingdon  
OX14 4SE, Oxon, UK  
Tel: +44 (0) 1235 828 200  
Fax: +44 (0) 1235 820 482

### Switzerland

Centro Nord-Sud 2E  
CH-6934 Bioggio-Lugano  
Tel: +41 (0) 91 604 55 22  
Fax: +41 (0) 91 605 17 85

### Deutschland

Bockenheimer Landstr. 17/19  
60325 Frankfurt/Main  
Tel: +49 (0) 69 779099  
Fax: +49 (0) 69 13376880

### United States

23591 El Toro Rd, Suite #167  
Lake Forest, CA 92630  
Tel: + 1 800 987 0985  
Fax: + 1 949 265 7703